



Win10 Driver Installation Guide

Index


1.Plug in the spectrometer.....	2
2.Disable driver signature enforcement	4
3.Intall Driver.....	8

1.Plug in the spectrometer

Right click to open [my computer](#) -> [computer management](#) -> [device manager](#) interface and plug in the spectrometer to the computer through USB. The following image will show up. It means the driver is not properly installed yet.

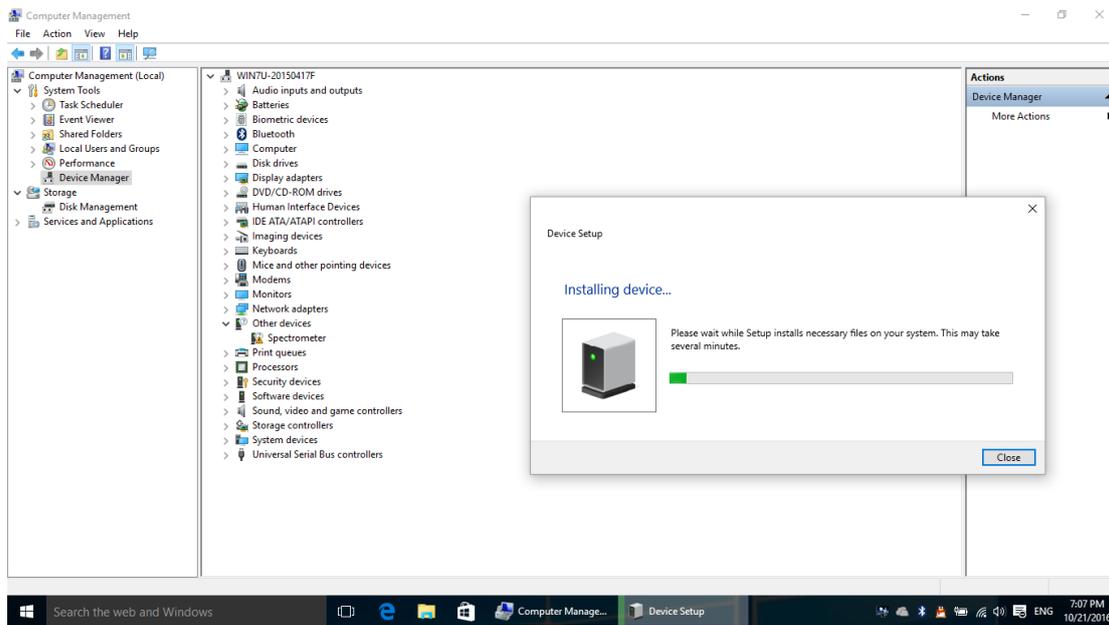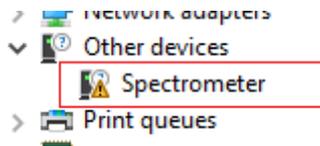

Since the driver does not contain digital signature, It needs to [disable driver signature enforcement](#) and then intall the driver properly.

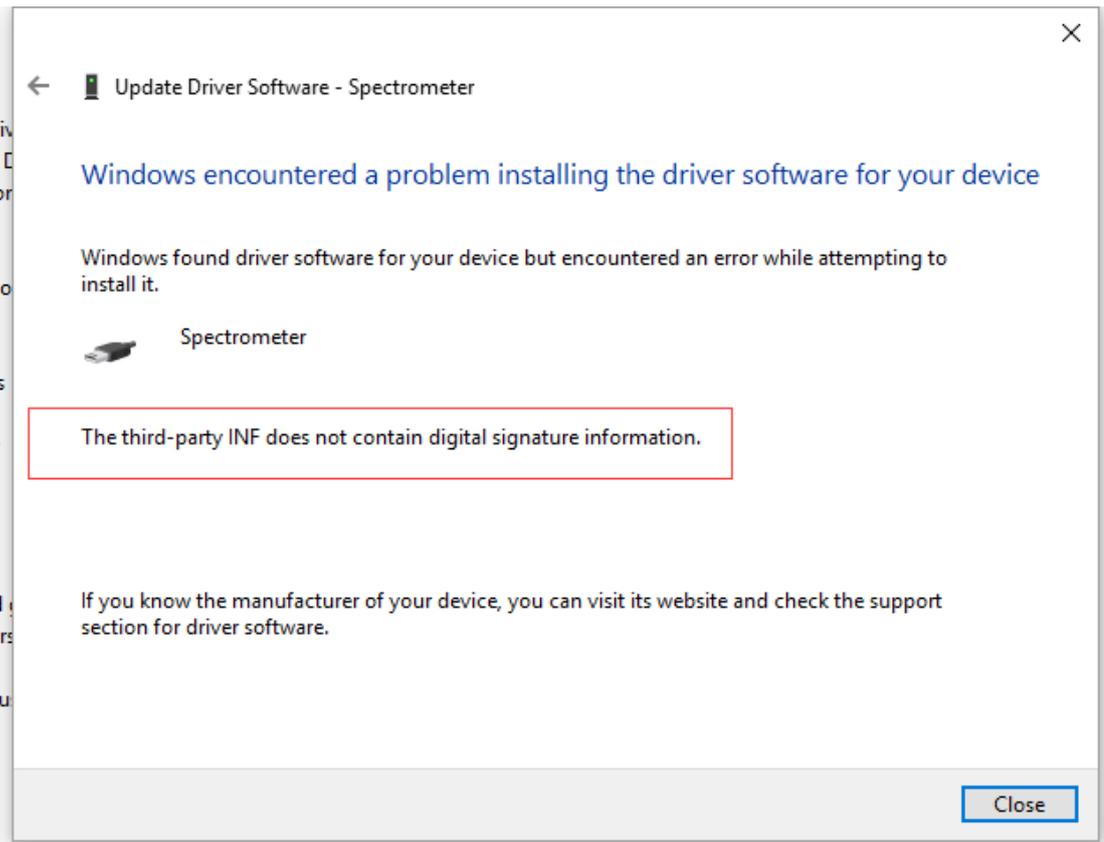


2.Disable driver signature enforcement


2.1 Click Windows start menu and enter the [Settings](#) interface




2.2 Enter "Update & security" under settings interface

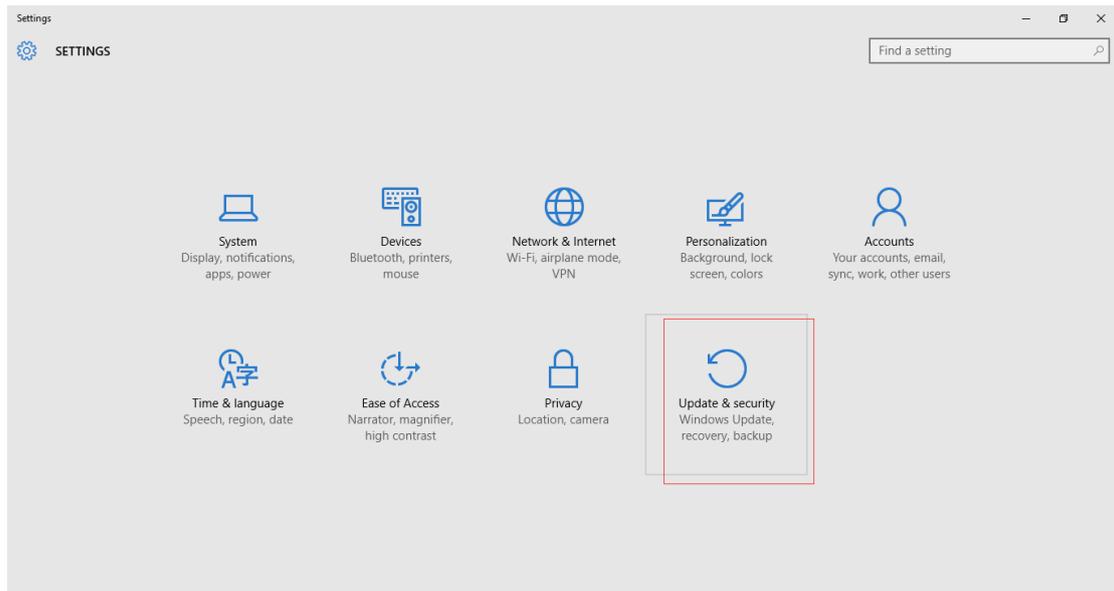


2.3 Find [Recovery](#) and click "[Restart now](#)" under the "[Advanced startup](#)" to restart the computer.

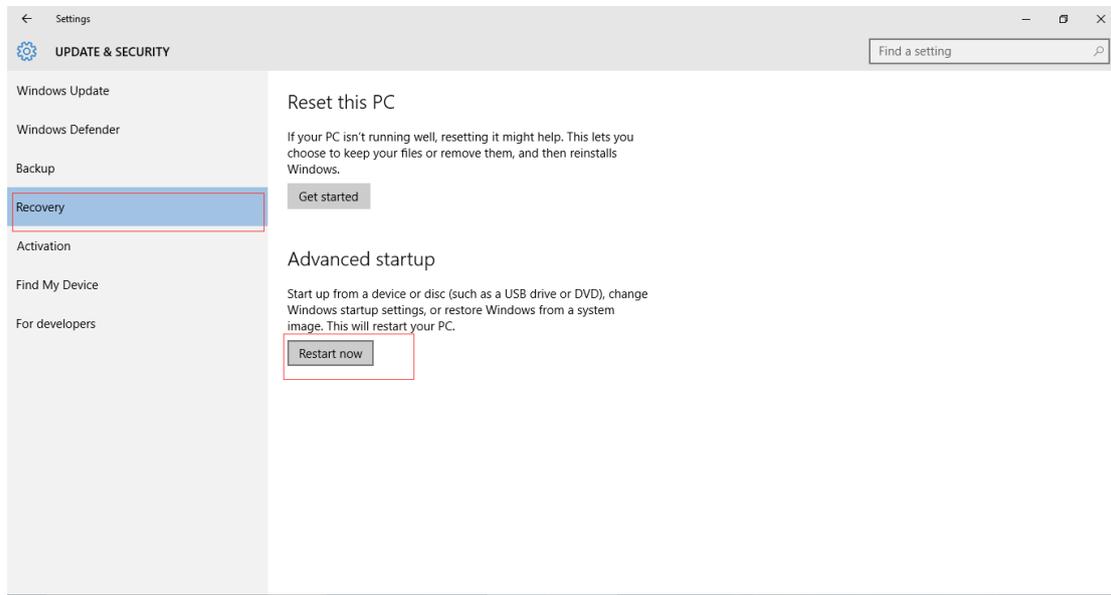


2.4 After the computer restart, choose "[Troubleshoot](#)".

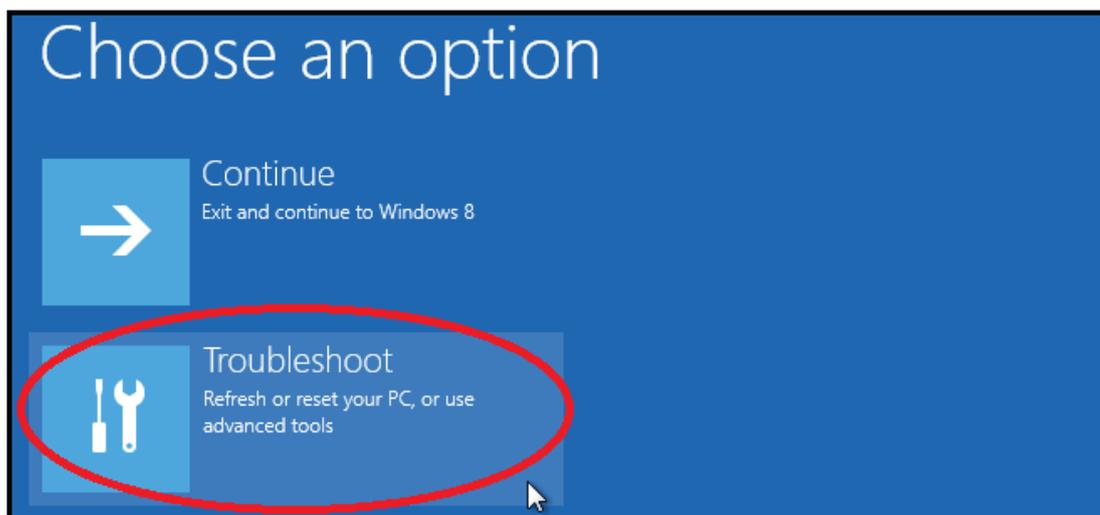


2.5 Select "[Advanced options](#)"

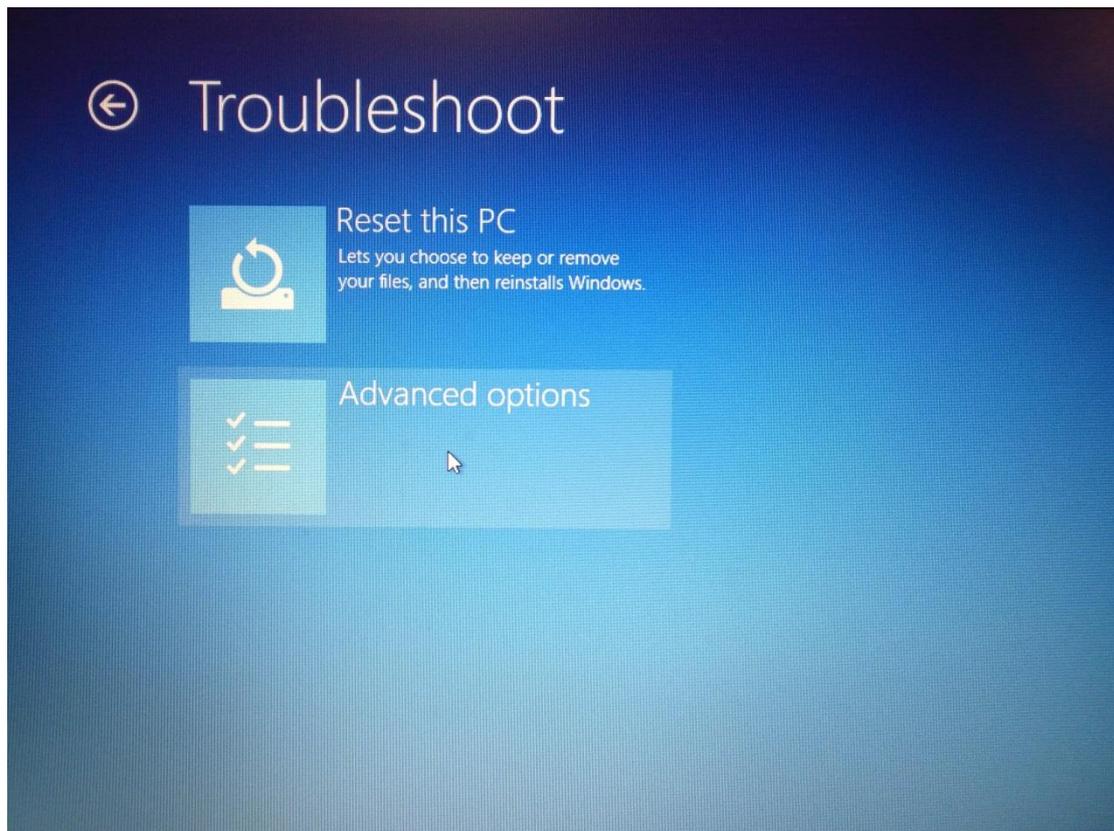


2.6 Select "[Startup settings](#)".

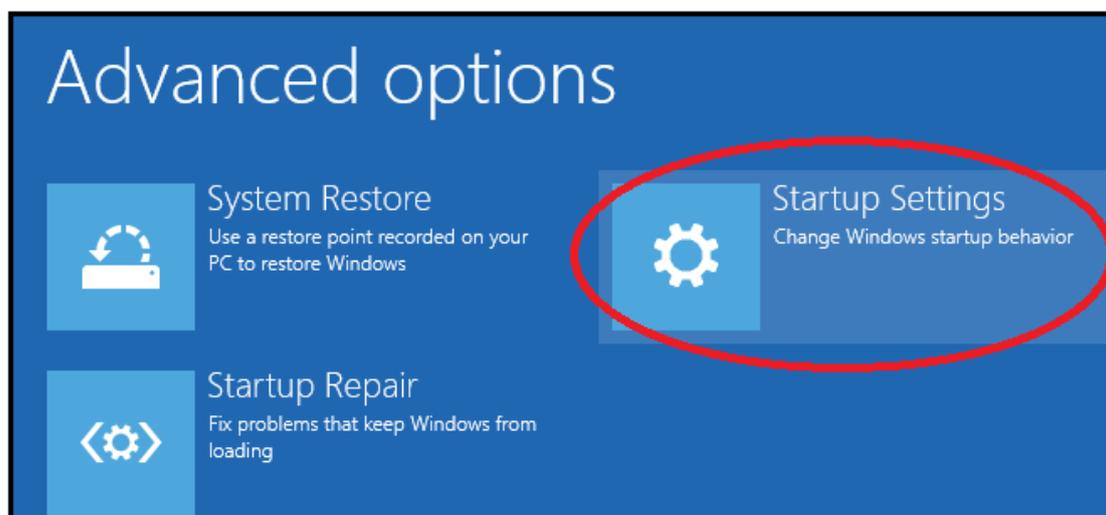

2.7 Click "Restart"


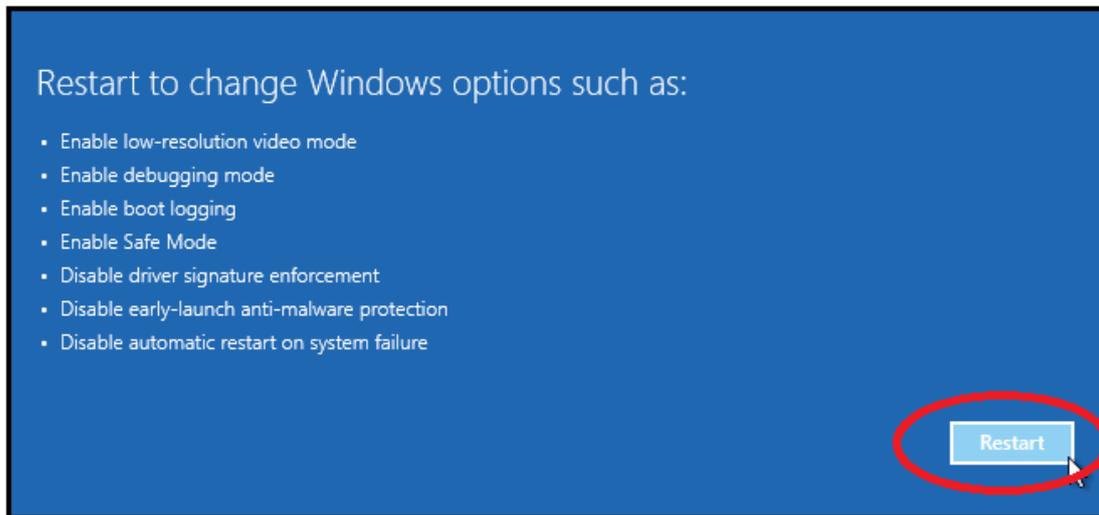



2.8 Select "7" to Disable Driver Signature Enforcement.

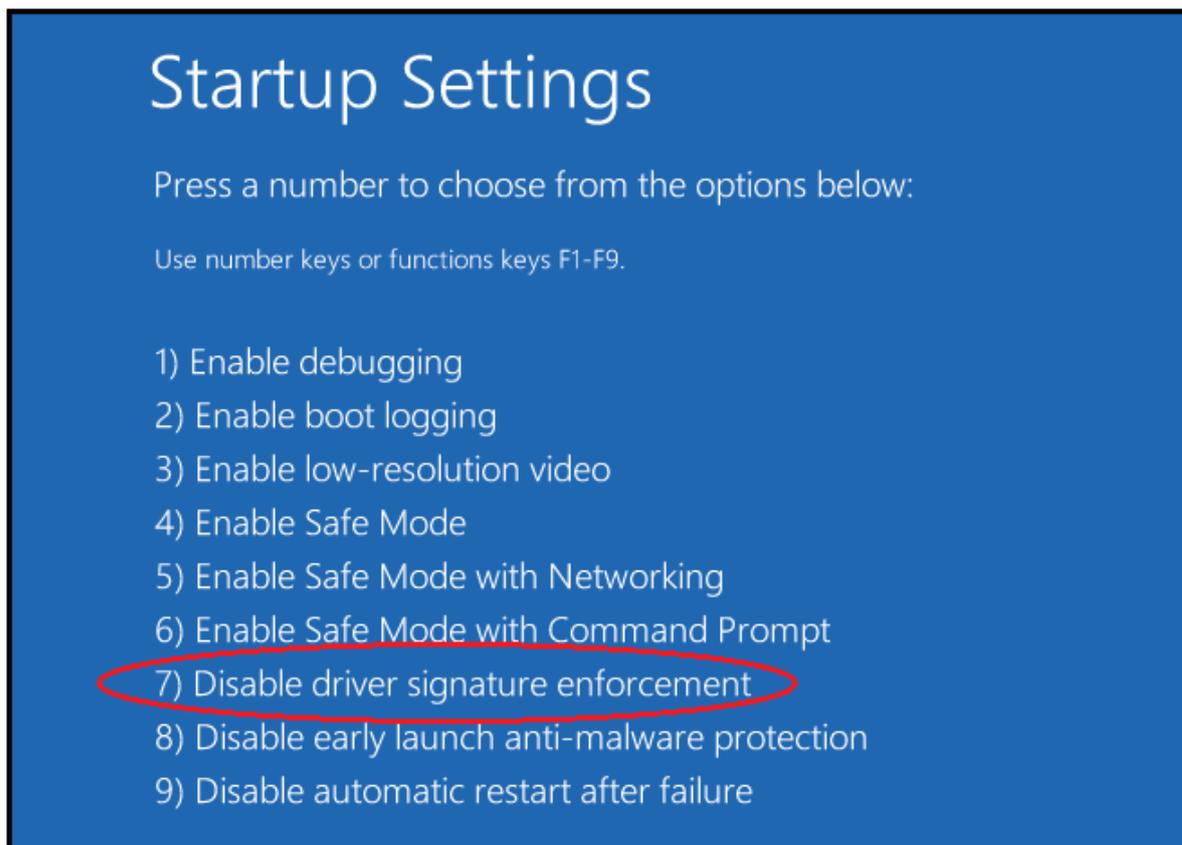

3.Intall Driver

Since the driver signature enforcement is disabled, we can intall the driver as follows:

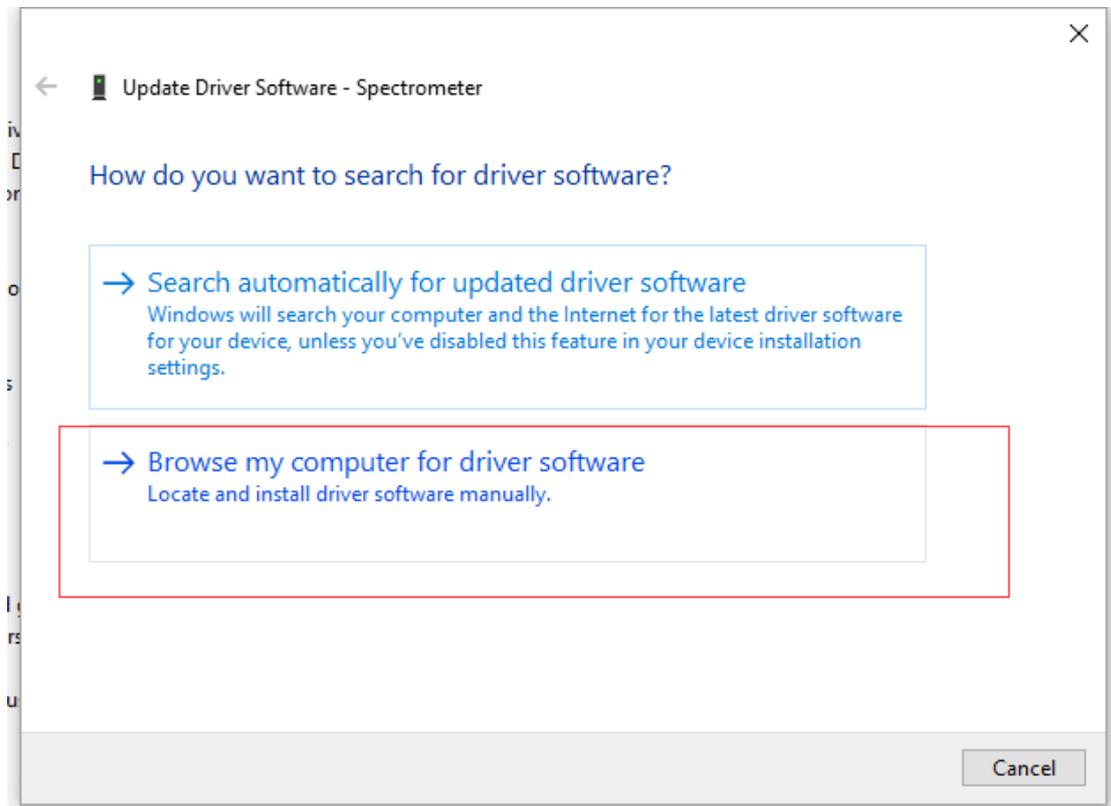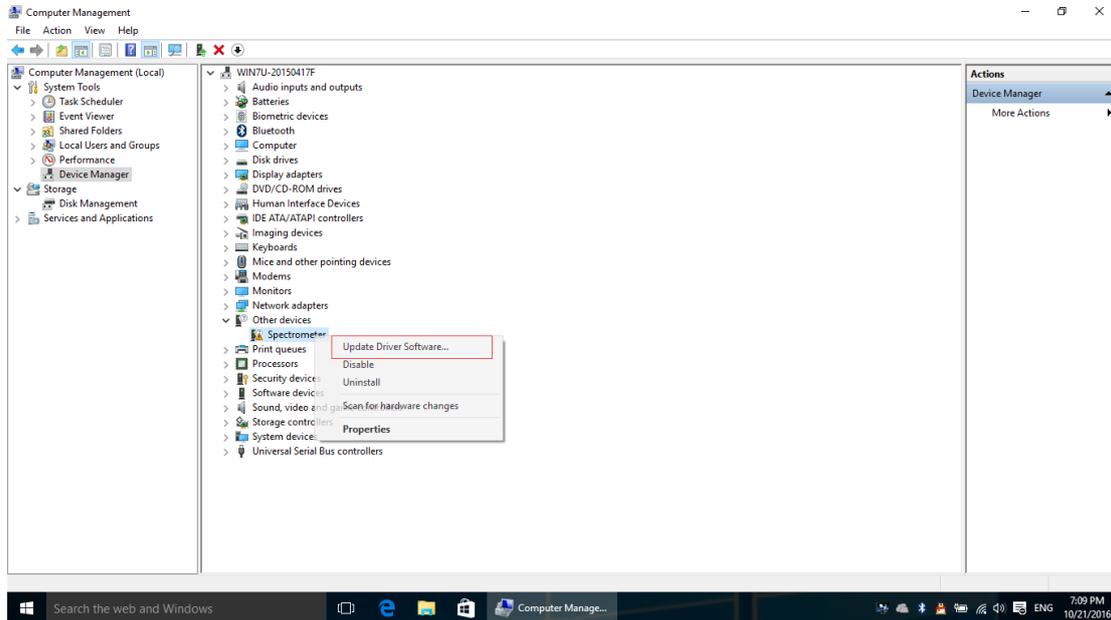

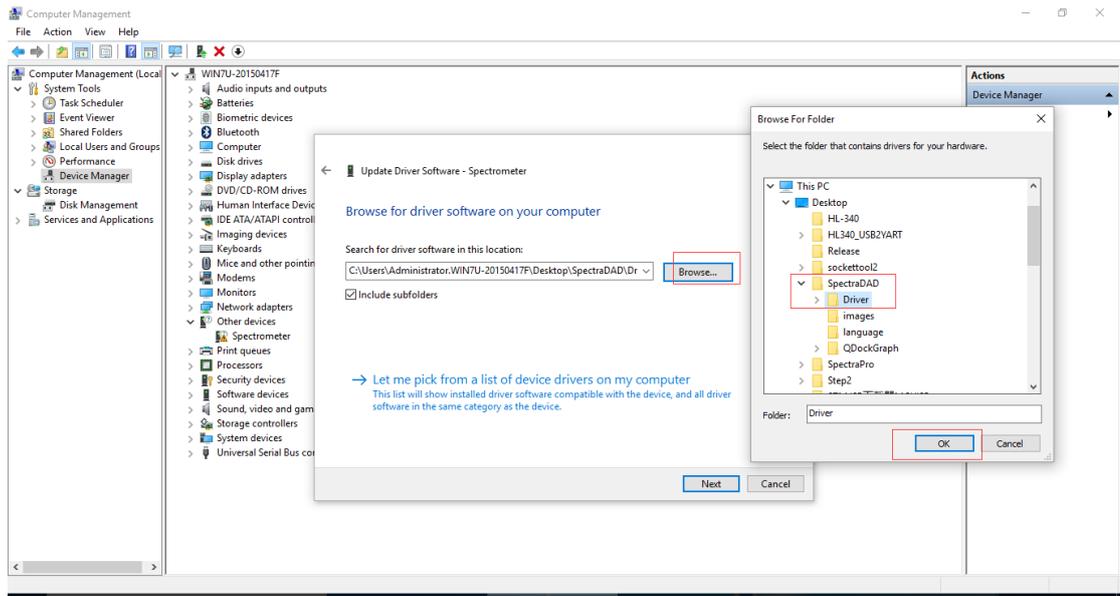

The driver files is located in the folder “**Driver**” after software installation.